

# AEROSIL® 200 Pharma – well established and ready for tomorrow's drug production

Continuous manufacturing is one of the greatest trends in the pharmaceutical industry, bringing advantages compared to the traditional batch production:

- Quality and cost position of the final medicine is improved
- Batch-to-batch variations are eliminated
- Intermediate storage becomes obsolete
- Potential future regulatory requirements are met.



As part of continuous manufacturing, continuous direct compression requires the integration of continuous powder feeding units, a continuous dry powder mixer and tablet press. Together, these production steps build the core equipment for a successful implementation of this manufacturing technique of growing importance.

Besides the clear advantages, continuous direct compression creates some challenges for the manufacturing process. Exact dosing of small amounts of Active Pharmaceutical Ingredients (APIs) and excipients is key to ensure a properly working continuous process. There is no downtime during the continuous run. The process must operate with low maintenance. Any issues, for example in powder flow, can impair the continuous production.



Gericke mixing equipment, with courtesy of Gericke GmbH

AEROSIL® Pharma is the industry standard glidant since decades. It ensures proper free flow within the tableting equipment and a homogeneous distribution of APIs. In addition, tablet properties such as mass variation and tablet hardness are improved by Evonik silica glidants. We at Evonik trust that continuous direct compression needs a glidant to be implemented successfully. This was demonstrated in an internal case study in cooperation with Gericke and L.B. Bohle.

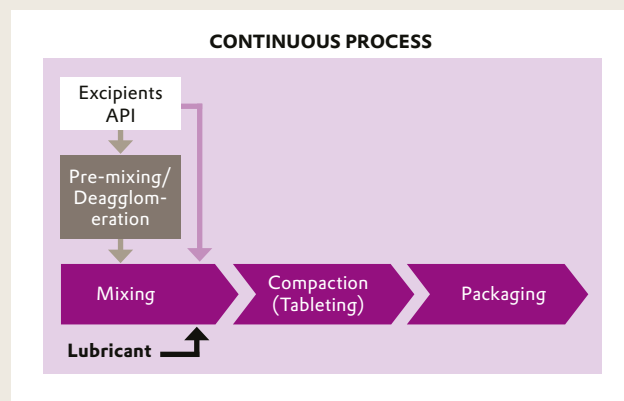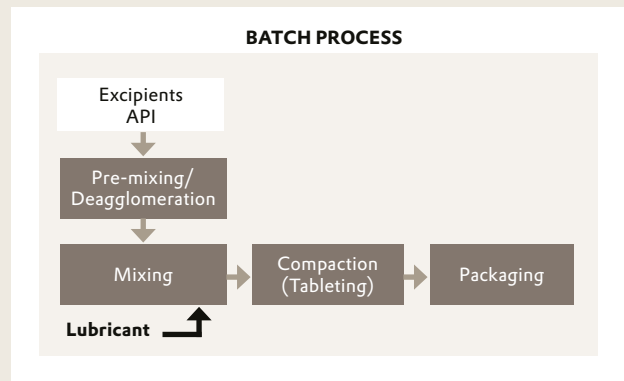

 Batch  Continuous

Direct Compression: Batch process vs. continuous process

The dosing of a glidant such as AEROSIL® 200 Pharma at very low feeding rates might be perceived as a challenge. With the proper feeding equipment, dosing of down to 100 g/h can be realized for the Evonik fumed silica.

The comparable low residence time of the materials in the continuous mixer has been shown to be sufficient to incorporate the AEROSIL® 200 Pharma into the system. The glidant is distributed on the surface of the host powder and can fulfill its function of decreasing the van der Waals forces, as well as, eliminating liquid bridges between particles in the mixture.

**INTERNAL FORMULATION CASE STUDY**

**Ingredient**


Microcrystalline Cellulose <sup>1</sup>	99.5 w.-%	99.0 w.-%
AEROSIL® 200 Pharma		0.5 w.-%
Magnesium stearate <sup>2</sup>	0.5 w.-%	0.5 w.-%

<sup>1</sup> Avicel PH 101, Sigma-Aldrich Chemie GmbH, Germany

<sup>2</sup> Magnesium Stearat, Caesar & Loretz GmbH, Germany

AEROSIL® 200 Pharma ensures the consistent powder flow from the mixer to the tableting machine. It reduces the retention of the powder in the tubing from the mixer and facilitates the filling of the compression chamber.

**TABLET MASS AND CRUSHING STRENGTH**

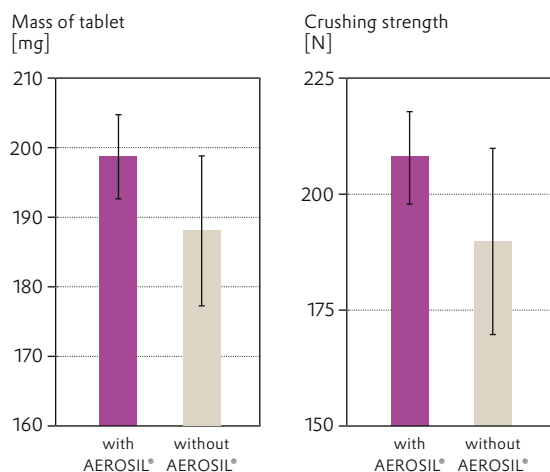

In the final tablets produced during the case study in the QbCon® System at L.B. Bohle, AEROSIL® 200 Pharma translates into higher tablet weights but strongly reduced tablet weight variation. In addition, the crushing strength is increased as well whereas its variance is decreased.



Dry Granulation line CM-Linie QbCon® DG at L.B. Bohle, with courtesy of L.B. Bohle Maschinen und Verfahren GmbH

Overall, AEROSIL® 200 Pharma is not only suitable for continuous direct compression, it is a central excipient for homogeneous and stable tablets produced in the continuous process, now and in future technologies.

This information and any recommendations, technical or otherwise, are presented in good faith and believed to be correct as of the date prepared. Recipients of this information and recommendations must make their own determination as to its suitability for their purposes. In no event shall Evonik assume liability for damages or losses of any kind or nature that result from the use of or reliance upon this information and recommendations. EVONIK EXPRESSLY DISCLAIMS ANY REPRESENTATIONS AND WARRANTIES OF ANY KIND, WHETHER EXPRESS OR IMPLIED, AS TO THE ACCURACY, COMPLETENESS, NON-INFRINGEMENT, MERCHANTABILITY AND/OR FITNESS FOR A PARTICULAR PURPOSE (EVEN IF EVONIK IS AWARE OF SUCH PURPOSE) WITH RESPECT TO ANY INFORMATION AND RECOMMENDATIONS PROVIDED. Reference to any trade names used by other companies is neither a recommendation nor an endorsement of the corresponding product, and does not imply that similar products could not be used. Evonik reserves the right to make any changes to the information and/or recommendations at any time, without prior or subsequent notice. AEROSIL® is a registered trademarks of Evonik Industries or its subsidiaries.

**Evonik Operations GmbH**  
Silica business line  
Rodenbacher Chaussee 4  
63457 Hanau  
Germany

Phone +49 6181 59-12532  
Fax +49 6181 59-712532  
ask-si@evonik.com  
[www.silica-specialist.com](http://www.silica-specialist.com)

**The Silica specialists at Evonik – Inside, to get it right.**